

- 
- 1) Write your name and record samples on the log sheet
  - 2) Log into the PC
  - 3) Open the WinUV/Scan software from the desktop (Cary WinUV>Scan)
  - 4) Once scan window is on the screen, turn on the power to the instrument and give it a minute or two to initialize (bottom lower left corner of the scan window shows status of the initialization)
  - 5) "Cary Online" should show up in top left corner of the Scan window when communication is successful. If "online" does not show up, close the scan window and re-open.
  - 6) Analyze samples
  - 7) When finished, turn off the lamps and leave the instrument on to allow the fan to cool it for about 5 minutes
  - 8) Shutdown the Cary WinUV software completely
  - 9) Turn off the power to the instrument