

HAAKE C10 Circulator Water Bath

Celeste Morin, x3756, Druckenmiller 243

November 14, 2017

Before Turning on the Water Bath:

- 1) Check to make sure the level of water in the bath is at an appropriate level.
 - a) Minimum: Up to 5cm below cover plate
 - b) Maximum: Up to 2cm below cover plate
 - c) Use DI water only to fill the bath
 - d) If water needs to be drained, scoop or siphon the water out of the bath.
This model does not have a drainage nozzle.
- 2) Check all hose connections to make sure they are all tight and none are cracking due to excess heat exposure.
- 3) The rotation speed of the pump motor can be changed with Switch (14).

